

Cedex Bio HT analyzer - MAINTENANCE SHEET

Serial No:

Year:



	Approximate system time (min)	Approximate operator time (min)	Jan	Feb	Mar	Apr	May	Jun	Jul	Aug	Sep	Oct	Nov	Dec
EVERY MONTH														
Clean Waste Box Fitting	-	1												
EVERY THREE MONTH **														
Clean external water reservoir and fluid waste reservoir	-	5												
Clean and lubricate Rotor	2	13												
Replace ventilation filters	-	5												
Replace external water reservoir filter	-	5												
EVERY SIX MONTH - Roche Service														
Clean Internal Water Reservoir	-	8												
Clean Wash Station	-	8												
Clean Instrument	-	16												
Roche Service	-	5												
WITHOUT INTERVAL - Status will be checked during Roche service visits														
Replace probe B / C	6	3												
Replace plunger tip dosage pipette B / C	5	5												
COUNTED INTERVAL - Usually done every six month during Roche service visits														
Replace Abs. Halogen lamp ***	21	5												
OPERATORS INITIALS														

** This actions are also part of the half year preventive maintenance. Hence, twice a year this actions are performed by the Roche Service.

*** 800h counter